



**Fast and Flexible Sample Preparation  
in Dioxin and PCB Analysis through Automation**

# Sample Preparation Done by Specialists

The automated DEXTech Plus sample preparation system for dioxin and PCB analysis cleans quickly and flexibly all kinds of extracts:

- food and feed samples, such as fish, meat, fish oil, eggs, vegetable oils and feed mixtures
- environmental samples, such as soil, sludge, sediments
- biological matrices, for example blood.

## „Greener Technology“

- Low solvent consumption
- Little waste (only glass, PTFE and filling materials)
- Reusable carbon column



## Advantages



### Fastest processing times

when processing from raw extracts (up to 5 g fat)



### Lowest costs



### Greatest flexibility

for various options see page 5



### Highest degree of automation





# Faster than Ever

3500 samples  
per year

Sample preparation with the novel DEXTech Plus 3-column-design is faster than ever. For example, a sample can be processed in only 35 min from raw extract; manual preparation is not necessary. Hence, only one system is required ( $\triangle$  one flow path) to process 3.500\* samples per year. Solvent consumption per run is very low.


Moreover, DEXTech Plus sets new standards in regards of flexibility, low cost per sample and effortless handling.

For high-throughput laboratories particularly attractive: Parallel use of 4 systems achieves a throughput of 14.000\* samples per year.



Set-up with four DEXTech Plus systems

\* Sample throughput with a changeover time of 3 minutes between samples, and a running time of 9 hours per day over 250 days per year

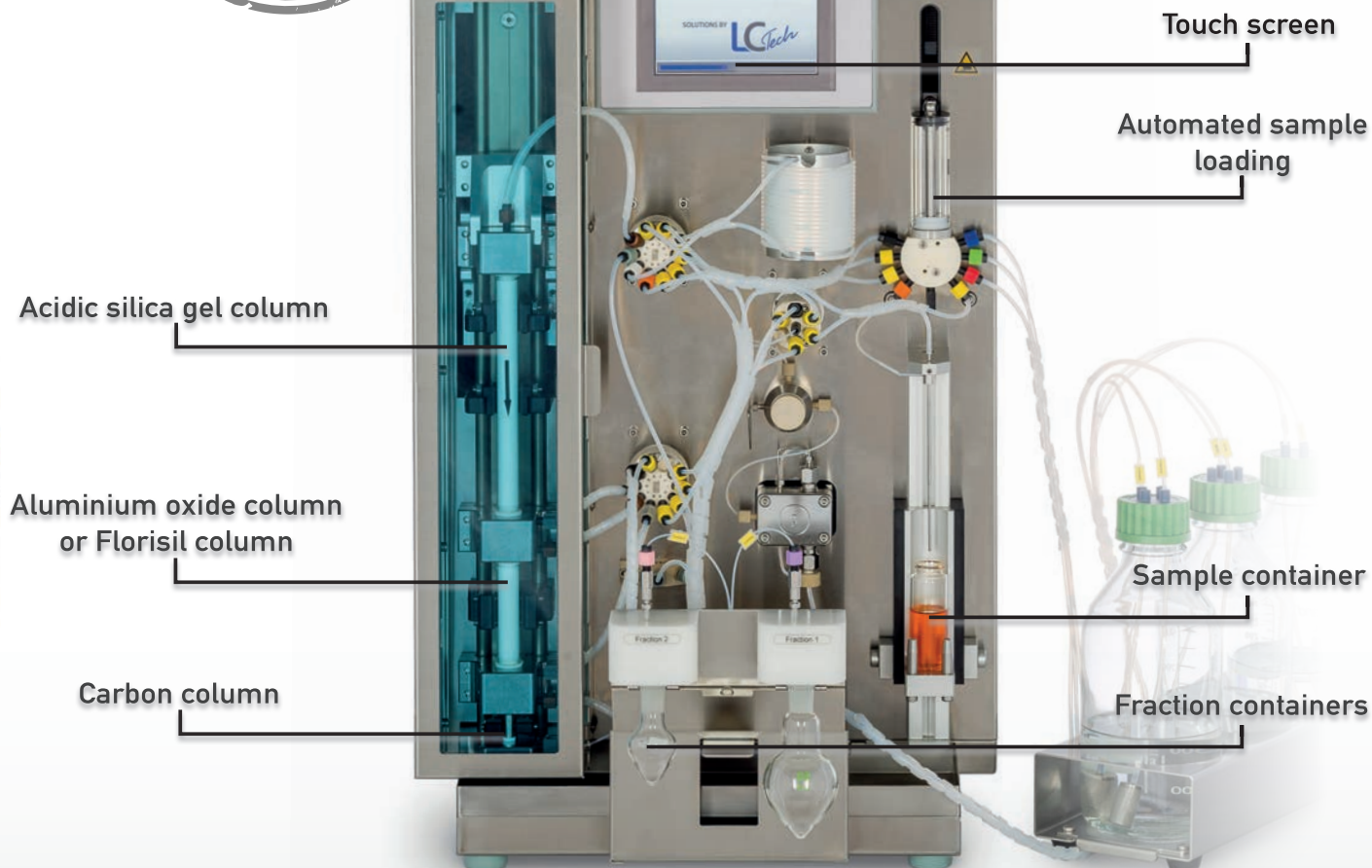

# Patented Locking System: Just Click It



## Low Cost 3-Column Set-up

### 1. Acidic silica gel column

- samples containing < 1.5 g fat – SMART column
  - samples containing > 1.5 - 5 g fat - Universal column
- Easy change between Universal and SMART column by fitting an adapter.

### 2. Aluminium oxide column or Florisil column;

depending on number and composition of target fractions.

### 3. Carbon column; reusable. This reduces the already low column cost even further.

Depending on the individual requirements, matrix or desired method, the user can equip the DEXTech Plus with different set-ups, simply insert the respective columns – „just click it“.

No screw connections, no tools, no other manual steps.  
From sample to sample – fast, flexible, effortless.

## Standard Method or Flexible Parameterisation

Evaluated standard methods, e.g. the default SMART method for 35-min processing, are stored in the DEXTech Plus. These methods can easily be accessed via the password-protected touch screen. The user may also freely parameterise a method and store it in the system. The stored methods can be opened and modified at any time and then be saved again under a different name.

Method Name	Default Smart	N°	1
<div> <div>back</div> <div>Conditioning</div> </div>			
	ml/min	min	n-Hexane
Conditioning 1	7.0	4.0	n-Hexane
Conditioning 2	7.0	2.0	n-Hexane
Conditioning 3	7.0	1.0	Toluene
Conditioning 4	7.0	1.0	DCM/n-Hexane

Screen for freely parametrizable method input




Easy, quick replacement and insertion of columns and adapter – „just click it“.



# Great Flexibility with Low Solvent Consumptions

## Aluminium Oxide or Florisil: The Choice Is Yours

By using different types of columns and the freely parametrizable method input, you have the opportunity to configure the DEXTech Plus system according to your requirements. The most common applications are shown below. Using the Florisil column you have the choice between two or three fractions, simply by changing the tubes.

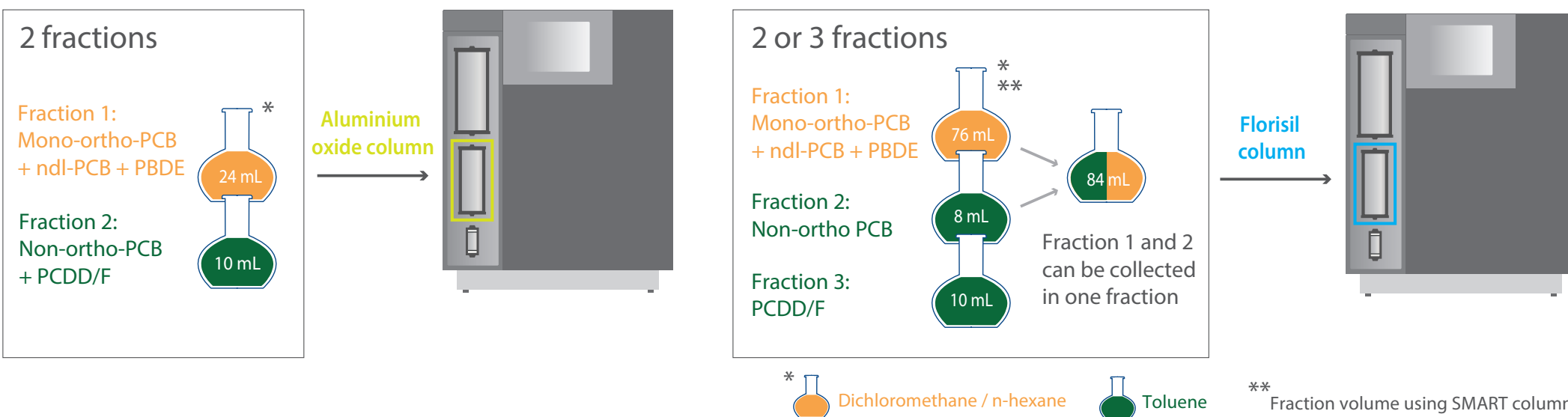

## Short Run Times, Low Solvent Consumption

Solvent	SMART column	Universal column
n-hexane	112 mL	203mL
DCM / n-hexane 1:1	24 mL	24 mL
Toluene	10mL	10 mL
<b>Total</b>	<b>146 mL</b>	<b>237 mL</b>

<b>Run time</b>	<b>34 min.</b>	<b>51 min.</b>
-----------------	----------------	----------------

Solvent	SMART column	Universal column
n-hexane	94 mL	178mL
DCM / n-hexane 1:1	24 mL	24 mL
Toluene	18mL	18 mL
<b>Total</b>	<b>136 mL</b>	<b>220 mL</b>

<b>Run time</b>	<b>44 min.</b>	<b>56 min.</b>
-----------------	----------------	----------------



# Automation with Safety



## Highest Degree of Automation

- Click columns into the column tower – “just click it”. The tower will be automatically locked by a patented process.
- Place sample vial into the DEXTech Plus. The sample will be automatically loaded via the syringe pump; however, at no time the sample will come in contact with the syringe pump itself. For quantitative transfer the sample container will be rinsed and the remains added into the sample loop.
- Select method and start process.

Throughout the entire process, the touch screen displays the processing status. An acoustic signal as well as an LED display indicates the end of processing.

The clean-up performed by the DEXTech Plus is compliant with the EPA method and European regulations.

## User Safety

- Leakage sensor below a removable drip tray
- Pressure sensor for pressure monitoring
- Overflow sensor for waste level control
- No mechanical movement when the door is open
- System operates in a low voltage range
- No fume hood necessary due to closed fraction flasks

Should one of the sensors send a signal, the system will stop processing. After elimination of the error, processing of the current sample will be continued.

## Report for Quality Control

As standard, DEXTech Plus is equipped with an USB interface to transfer the in the software generated processing protocols to a PC. All sample related data are consequently stored for documentation.




*USB interface for documentation*



*Automated sample loading*




*Rinsing of sample vial for quantitative transfer*

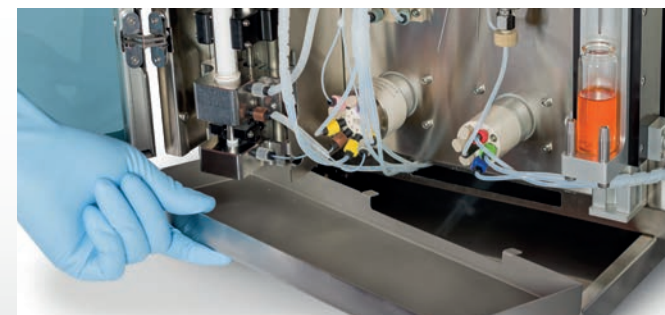

*Easy to remove and clean drip tray above the leakage sensor*




# Reliable and Reproducible Results

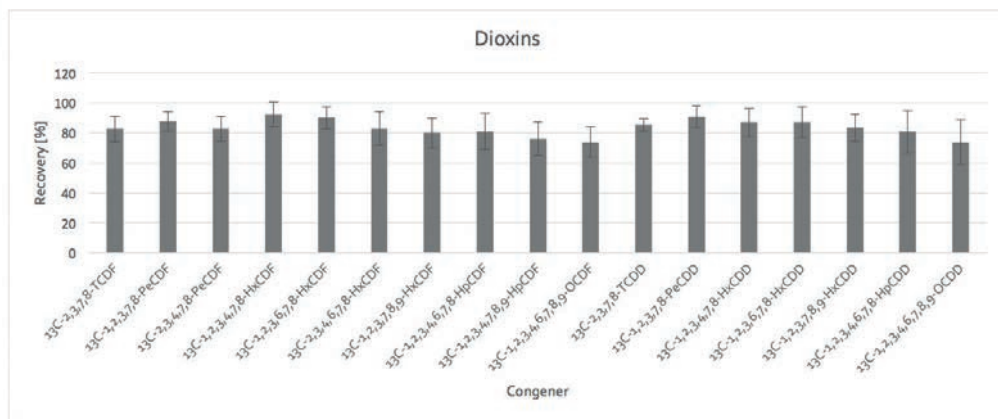

## Results Dioxin

Sources: Chemical and Veterinary Investigation Office Münsterland-Emscher-Lippe (governmental laboratory); LCTech Laboratory

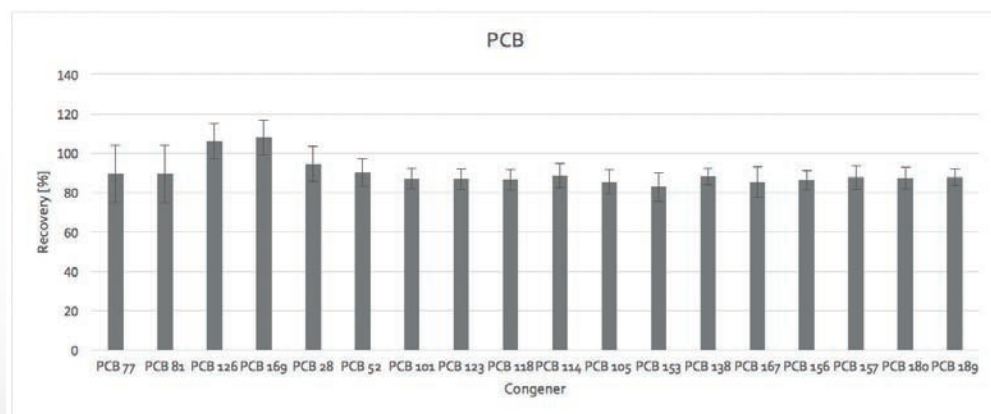


## Results PCB

Sources: Chemical and Veterinary Investigation Office Münsterland-Emscher-Lippe (governmental laboratory); LCTech Laboratory

## No Cross-Contamination

Cross-contamination is reliably prevented.

- Continuous rinsing of all parts that have come in contact with the sample
- Exchangeable columns
- No contact between sample and syringe pump

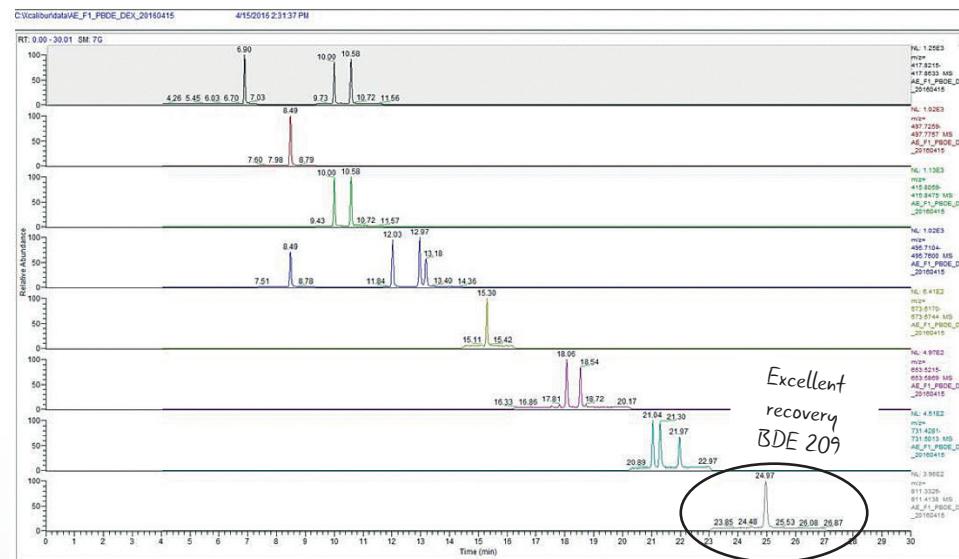

Clear signals, sharp separation, low background noise and excellent recoveries

Sources: Bavarian Health and Food Safety Authority (LGL-Oberschleißheim, governmental laboratory); LCTech Laboratory

## We are proud to compare:



Lowest  
Cost



Fastest  
Processing Time



No Cross-  
Contamination



Highest Degree  
of Automation



Greatest  
Flexibility



Ultimate  
User Safety



The information contained in this brochure is based on our current knowledge and has been carefully checked. However, since we continually work on the further development of our products, please accept texts, pictures and numbers on these pages as non-binding and exemplary only.

SOLUTIONS BY **LC***Tech*

LCTech GmbH  
Daimlerstr. 4, 84419 Obertaufkirchen, Germany  
Tel: +49 8082 2717-0 Fax: +49 8082 2717-100  
dioxins@LCTech.de [www.LCTech-online.com](http://www.LCTech-online.com)

